

Motion Detector 2




(Order Code MD-BTD)

The Motion Detector is used to collect position, velocity and acceleration data of moving objects. Students can study a variety of motions with the Motion Detector, including

- Walking toward and away from the Motion Detector.
- Dynamics carts moving on track.
- Objects in simple harmonic motion, such as a mass hanging on a spring.
- Pendulum motions.
- Objects dropped or tossed upward.
- A bouncing object.




This is motion data collected with Logger Pro® and a computer.

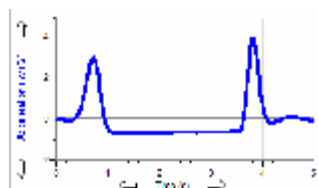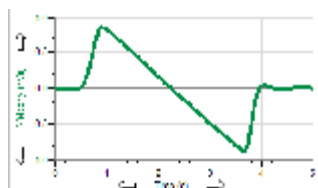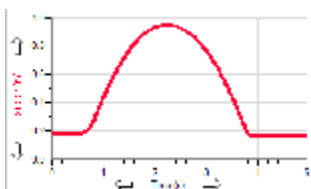


Ball Toss with Logger Pro


Here are some sample calculator data collected using a Motion Detector:

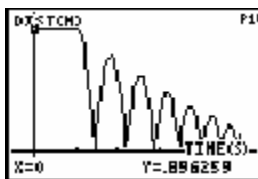

Ball Bouncing

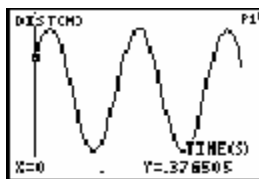

Weight on Spring

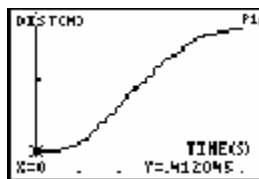

Student Walking

NOTE: This product is to be used for educational purposes only. It is not appropriate for industrial, medical, research, or commercial applications.

The Motion Detector is compatible with the following data-collection interfaces:

- Vernier LabPro®
- Texas Instruments CBL 2™
- Texas Instruments CBL™
- Universal Lab Interface (requires separate cable, order code ULI-MDC)


What is Included with the Motion Detector?

The Motion Detector ships with the detector itself and a cable that connects the Motion Detector to the LabPro, CBL 2 or original CBL interface.



How the Motion Detector Works

This Motion Detector emits short bursts of ultrasonic sound waves from the gold foil of the transducer. These waves fill a cone-shaped area about 15 to 20° off the axis of the centerline of the beam. The Motion Detector then “listens” for the echo of these ultrasonic waves returning to it. The equipment measures how long it takes for the ultrasonic waves to make the trip from the Motion Detector to an object and back. Using this time and the speed of sound in air, the distance to the nearest object is determined.

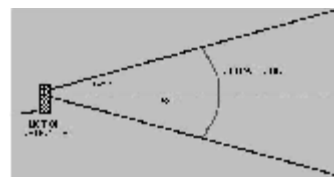

Note that the Motion Detector will report the distance to the closest object that produces a sufficiently strong echo. The Motion Detector can pick up objects such as chairs and tables in the cone of ultrasound.

The sensitivity of the echo detection circuitry automatically increases, in steps, every few milliseconds as the ultrasound travels out and back. This is to allow for echoes being weaker from distant objects.

Features of the Motion Detector

- The Motion Detector is capable of measuring objects as close as 0.15 m and as far away as 6 m. The short minimum target distance (new to this version of the Motion Detector) allows objects to get close to the detector, which reduces stray reflections.
- The Motion Detector has a pivoting head, which helps you aim the sensor accurately. For example, if you wanted to measure the motion of a small toy car on an inclined plane, you can lay the Motion Detector on its back and pivot the Motion Detector head so that it is perpendicular to the plane.
- The Motion Detector has a Sensitivity Switch, which is located under the pivoting Motion Detector head. To access it, simply rotate the detector head away from the detector body.



- Rubber feet on the Motion Detector provide stability. Rubber feet are located on the back of the Motion Detector, which provides stability when the Motion Detector is placed on its back. Rubber feet on the bottom provide stability when the Motion Detector is standing up.
- A standard camera screw mount can be found on the back of the Motion Detector. This screw mount allows you to use the Motion Detector Clamp (sold separately, order code MD-CLAMP, \$6), which is a universal-mounting clamp, or to use it with typical tripod-mounting hardware.



Camera Screw Mount



Motion Detector Clamp (sold separately)

Connecting the Motion Detector to an Interface


To use the Motion Detector with LabPro, CBL 2 or original CBL, the Motion Detector is connected to the interface with the included cable. Plug the cable into the DIG/SONIC port on the right side of the Motion Detector. Plug the other end of the cable into the interface. To use the Motion Detector with the Universal Lab Interface, plug the cable into Port 2 of the ULI.



Using the Sensitivity Switch

Slide the Sensitivity Switch to the right to set the switch to the “normal” setting. This setting is best used for experiments such as studying the motion of a person walking back and forth in front of the Motion Detector, a ball being tossed in the air, pendulum motion, and any other motion involving relatively large distances or with objects that are poor reflectors, e.g., coffee filters.

The other sensitivity setting, which we call “Track”, works well when studying motion of carts on tracks like the Vernier Dynamics System, or motions in which you want to eliminate stray reflections from object near to the sensor beam.

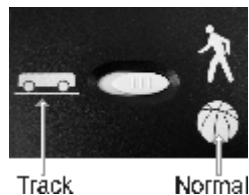

Using the Motion Detector with a Computer

This sensor can be used with a Macintosh® or PC computer and the Vernier LabPro. Here is the general procedure for using the Motion Detector with a computer:

1. Connect the Motion Detector to the DIG/SONIC port on the interface.
2. Start the *Logger Pro* software on the computer.
3. You may alternately open an experiment file for the Motion Detector, and you are ready to collect data.

Using the Motion Detector with TI Graphing Calculators

This sensor can be used with a TI graphing calculator and any of the following lab interfaces: LabPro, CBL 2™, or CBL™. Here is the general procedure to follow when using the Motion Detector with a graphing calculator:

1. Make sure that a data-collection program is installed on your calculator. If you are using LabPro or CBL 2, use the DataMate program. If you are using the original CBL, use the Physics program.
2. Use the calculator-to-calculator link cable to connect the interface to the TI graphing calculator using the I/O ports located on each unit. Be sure to push both plugs in firmly.
3. Connect the Motion Detector to any of the sonic ports on the interface.
4. Start the data-collection program, and you are ready to collect data.

Using the Motion Detector with Palm OS® Handhelds

This sensor can be used with a Palm OS handheld and the LabPro.

1. Connect the Palm OS handheld, LabPro, and the Motion Detector.
2. Start Data Pro.
3. Tap New, or choose New from the Data Pro menu. Tap New again. The Motion Detector will be identified automatically.
4. You are now ready to collect data.

Using the Motion Detector with Other Sensors

The Motion Detector can be used with one or more other sensors connected to the interface. Here are some examples of how the Motion Detector can be used for experiments with another sensor:

- With a force sensor to study the relationship between force and motion
- With a force sensor to study collisions and impulse
- With a force sensor to study simple harmonic motion
- With a light sensor to study the inverse square law
- With a magnetic field sensor to study how magnetic field varies with position
- With a second Motion Detector (LabPro only)

Tips on Getting Good Results with the Motion Detector¹

The most frequently reported problems with a Motion Detector are (1) that the Motion Detector does not work beyond a certain distance or (2) that the graph is very noisy. There are a number of ways to troubleshoot these situations, and they are described below.

1. See if the Sensitivity Switch makes a difference. Simply set the Sensitivity Switch to the other position and retry the experiment. This change may solve the problem.
2. The Motion Detector does not work beyond a certain distance, e.g., it does not detect anything beyond 1.2 m. Here are some things to check if you have this problem:
 - Check for movable objects (textbooks, ring stands, etc.) in the cone of the ultrasound. If possible, move these objects out of the measurement cone. It may not take a very large object to cause problems.
 - Check for a stationary object (chair, table, etc.) in the cone of the ultrasound. This object may be detected when you are trying to study an object further away. It may not take a very large object to cause problems. If you have trouble with a stationary object causing unwanted echoes, try setting the equipment up so that the objects are not in the coned or placing a cloth over the object. This minimizes the ultrasound reflection.
 - Also note that the cone of ultrasound extends downward from the center line. This can cause problems if you are using the Motion Detector on a hard, horizontal surface. In these cases, try pivoting the head of the Motion Detector to aim it slightly upward.
3. Noisy or erratic data may have a number of causes. Here are some tips.
 - Sometimes other sound sources can cause problems. If there is another source of ultrasonic waves in the same frequency range, this will cause erroneous readings. Examples include motors and fans, air track blowers, the sound made by the air exiting the holes on an air track, etc. Try to eliminate these sources of noise. If you are using an air track, try changing the air flow volume.
 - Make sure that the Motion Detector is not placed close to a computer or computer monitor.



- If the room in which the Motion Detector is being used has a lot of hard, sound-reflecting surfaces, you can get strange effects caused by the ultrasound bouncing around the room. Standing waves can be set up between the Motion Detector and a sound reflector. Try placing a cloth horizontally just in front of and below the Motion Detector. This sometimes helps eliminate ultrasound that is “skipping” into the Motion Detector.
- Try changing the data collection rate. Sometimes Motion Detectors work better at one data rate than another. Rates above 30 Hz do not work well in acoustically live rooms.
- If you are studying people moving, have them hold a large, flat object (e.g., a large book or a pizza box) as a reflector. If you have an irregular reflecting surface, sometimes the waves will be reflected back to the transducer, and sometimes not. The results will seem erratic.

Specifications

Range: 0.15 to 6 m

Resolution: 1 mm

Support Products

Motion Detector Clamp (order code MD-CLAMP, \$6)

The Motion Detector Clamp attaches to the back of the Motion Detector, and then the clamp can be attached to a variety of objects such as table tops, ring stands, etc. The clamp adds versatility



Vernier Dynamics System (order code VDS)


The Vernier Dynamics System consists of a versatile track and two carts with ultra low friction bearings. This system facilitates performing kinematics and dynamics experiments. The Motion Detector attaches to the dynamics track in such a way that the entire track can be used for motion studies.



¹ An excellent discussion of motion detector theory and operation can be found in “Physics and Technical Characteristics of Ultrasonic Sonar Systems”, Dan MacIsaac and Ari Hamalainen, *The Physics Teacher* **40**, 39 – 46 (January 2002).

Warranty

Vernier warrants this product to be free from defects in materials and workmanship for a period of five years from the date of shipment to the customer. This warranty does not cover damage to the product caused by abuse or improper use.

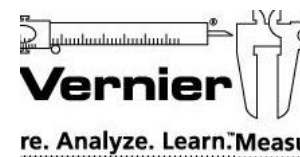

Vernier Software & Technology

13979 S.W. Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440
info@vernier.com • www.vernier.com

Rev. 10/26/05

Logger *Pro*, Vernier LabPro and other marks shown are our registered trademarks in the United States.
CBL 2 and CBL are trademarks of Texas Instruments.

All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.